



## BugDorm-4E3030D Specimen Handling Cage [BD4E3030D]

MegaView Science Co., Ltd.

<https://shop.bugdorm.com> · <https://megaview.com.tw>



### Package Contents

- x1 Fabric Cage Body
- x12 Fiberglass Rods
- x4 Three-way Joints
- x4 Webbed Three-way Joints

### Assembly

1. Thread one rod into each of the 12 sleeves of the cage.
2. Connect rods with the three-way joints. The webbed joints are designed to secure the hard floor (made of acrylic, sold separately). If the solid bottom is not required, all three-way joints are interchangeable and can be used at any corner to connect the rods.


### Note

- \* Operating Temperature: -10 to 60°C
- \* The soft plastic floor is not waterproof.
- \* Cage fabric is not UV treated. Avoid long-term exposure to direct sunlight.

**CLEANING** HAND WASH: Submerge and soak the cage for 20-30 minutes in the mix of detergent and water, agitate the cage gently, and then rinse the cage to remove the detergent.

MACHINE WASH: Use the side-loading laundry machine set on a washing cycle of low agitation and low-speed spin.

**CAUTION**: Wash cage body only. Wash at room temperature. Use laundry bag when machine wash. Recommend applying neutral detergents. Apply bleach or stain remover with care, as those chemicals weaken the cage fabric. Dry the cage in the shade.

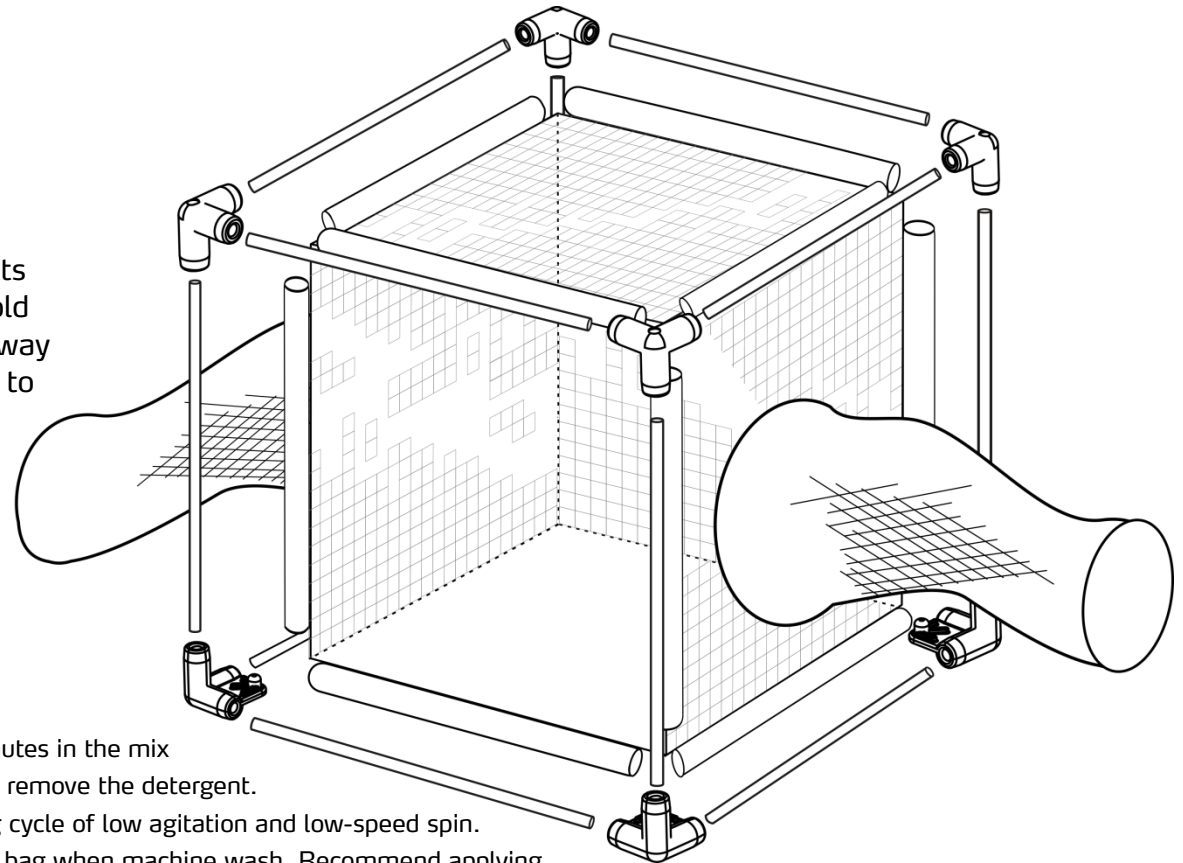

Have any questions? Please e-mail us: [bugdorms@megaview.com.tw](mailto:bugdorms@megaview.com.tw) or [bugdorms@gmail.com](mailto:bugdorms@gmail.com)

R20240710