



## BugDorm-4S4590DH Specimen Handling Cage [BD4S4590DH]

MegaView Science Co., Ltd.

<https://shop.bugdorm.com> · <https://megaview.com.tw>



### Package Contents

- x1 Fabric Cage Body
- x16 Fiberglass Rods
- x4 Two-way Joints
- x4 Three-way Joints
- x4 Webbed Three-way Joints

### Assembly

1. Use four two-way joints to connect eight rods into four long rods.
2. Thread one long rod from Step 1 into the two sleeves on each of the four long sides of the cage.
3. Thread the remaining eight rods into the short sleeves of the cage.
4. Connect rods with three-way joints. The webbed joints are designed to secure the hard floor for other cage models. They are interchangeable with ordinary joints and can be used at any corner to connect the rods.


### Note

- \* Operating Temperature: -10 to 60°C
- \* The soft plastic floor is not waterproof.
- \* Cage fabric is not UV treated. Avoid long-term exposure to direct sunlight.

**CLEANING** HAND WASH: Submerge and soak the cage for 20-30 minutes in the mix of detergent and water, agitate the cage gently, and then rinse the cage to remove the detergent.

MACHINE WASH: Use the side-loading laundry machine set on a washing cycle of low agitation and low-speed spin.

**CAUTION**: Wash cage body only. Wash at room temperature. Use laundry bag when machine wash. Recommend applying neutral detergents. Apply bleach or stain remover with care, as those chemicals weaken the cage fabric. Dry the cage in the shade.

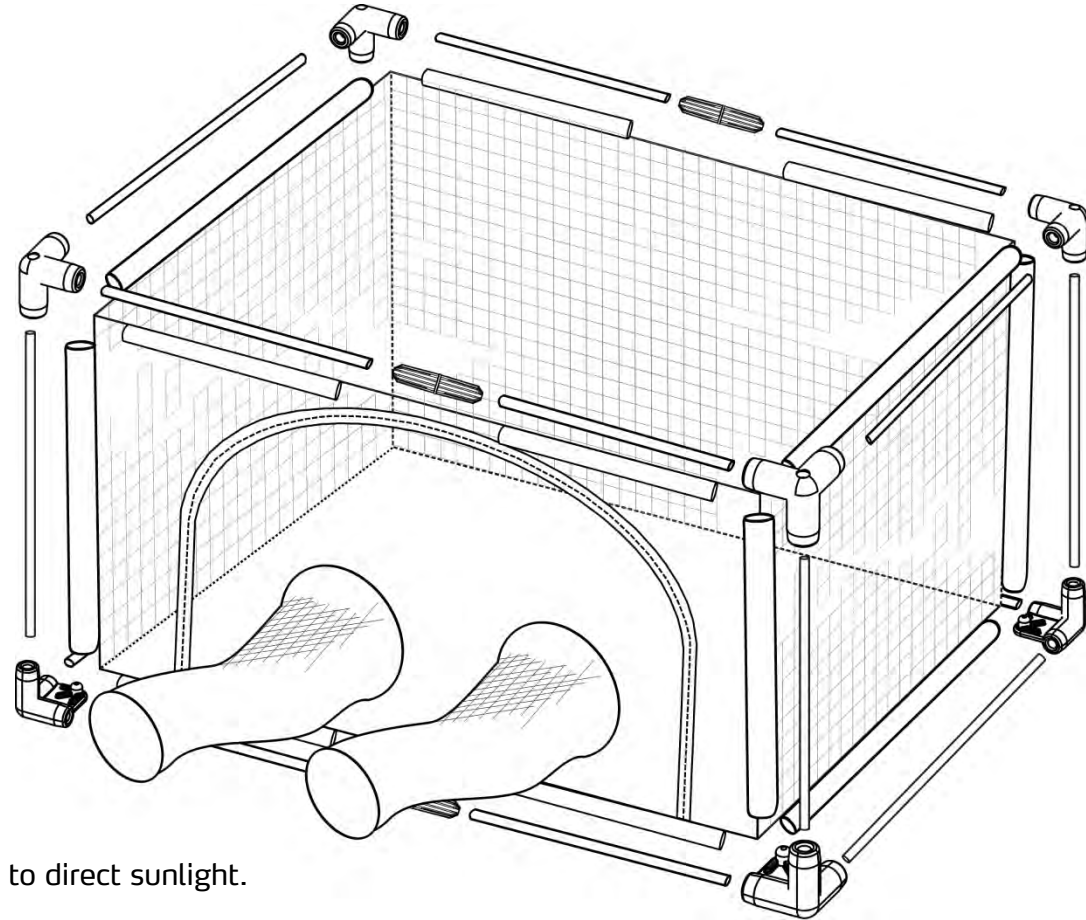

Have any questions? Please e-mail us: [bugdorms@megaview.com.tw](mailto:bugdorms@megaview.com.tw) or [bugdorms@gmail.com](mailto:bugdorms@gmail.com)

R20240710