



BugDorm Clip Cage ☐ BDC1540 ☐ BDC1540H ☐ BDC1560 ☐ BDC1560H

MegaView Science Co., Ltd.

<https://shop.bugdorm.com> · <https://megaview.com.tw>



Package Contents

Clip Cage [BDC1540] (pack of 12, ID*: 2.5cm)

x24 Foam Rings
x30 Mesh Circles (6 spares)
x48 Stainless Pins



Clip Cage with Access Hole [BDC1540H] (pack of 12, ID*: 2.5cm)

x12 Foam Rings
x12 Foam Rings with Access Hole
x30 Mesh Circles (6 spares)
x48 Stainless Pins



Clip Cage [BDC1560] (pack of 12, ID*: 4cm)

x24 Foam Rings
x30 Mesh Circles (6 spares)
x72 Stainless Pins



Clip Cage with Access Hole [BDC1560H] (pack of 12, ID*: 4cm)

x12 Foam Rings
x12 Foam Rings with Access Hole
x30 Mesh Circles (6 spares)
x72 Stainless Pins




*ID = Internal Diameter

Assembly



1. Peel off protective paper.

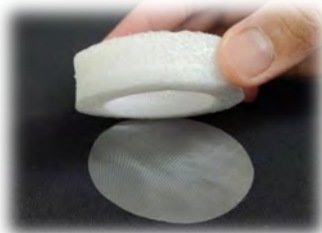

2. Press foam over mesh circle.



3. Trim off excess mesh.
Repeat Step 1 to 3
with another foam ring.



4. Sandwich leaf between
two foams.



5. Clamp foam rings
together with pins



6. A complete clip cage.



7. Pull off plug to access cage interior.
Available for H-type clip cages only.

Have any questions?

Please e-mail us: bugdorms@megaview.com.tw or bugdorms@gmail.com