

# Pop-up Berlese Funnel [BT5003]

MegaView Science Co., Ltd.

<https://shop.bugdorm.com> · <https://megaview.com.tw>





## Package Contents

- x1 Fabric Funnel Body
- x1 Centrifuge Tube (50 ml)
- x1 Tube Connector (2 parts)
- x1 Carrying Bag


## Assembly

The pop-up Berlese funnel is assembled out of the box. If you need to attach the fabric funnel to the tube connector again, please follow the steps below.

1. Thread the narrow end of the black funnel cloth through the upper piece of the connector.
2. Reverse the black cloth to surround the short cylinder of the upper piece of the connector.
3. Push and snap the upper piece of the connector in Step 2 into the bottom piece of the connector.
4. Push the centrifuge tube onto the bottom piece of the connector to complete the assembly.
5. If the connector pieces don't securely lock, use rubber bands, screws, or zip ties (Pic. A & B, NOT included) to tighten them up together.

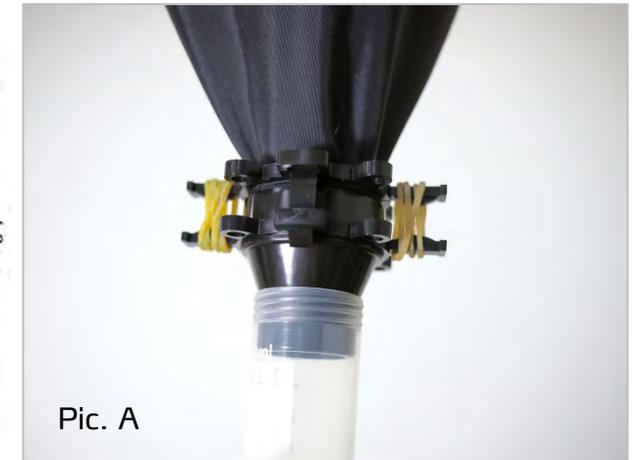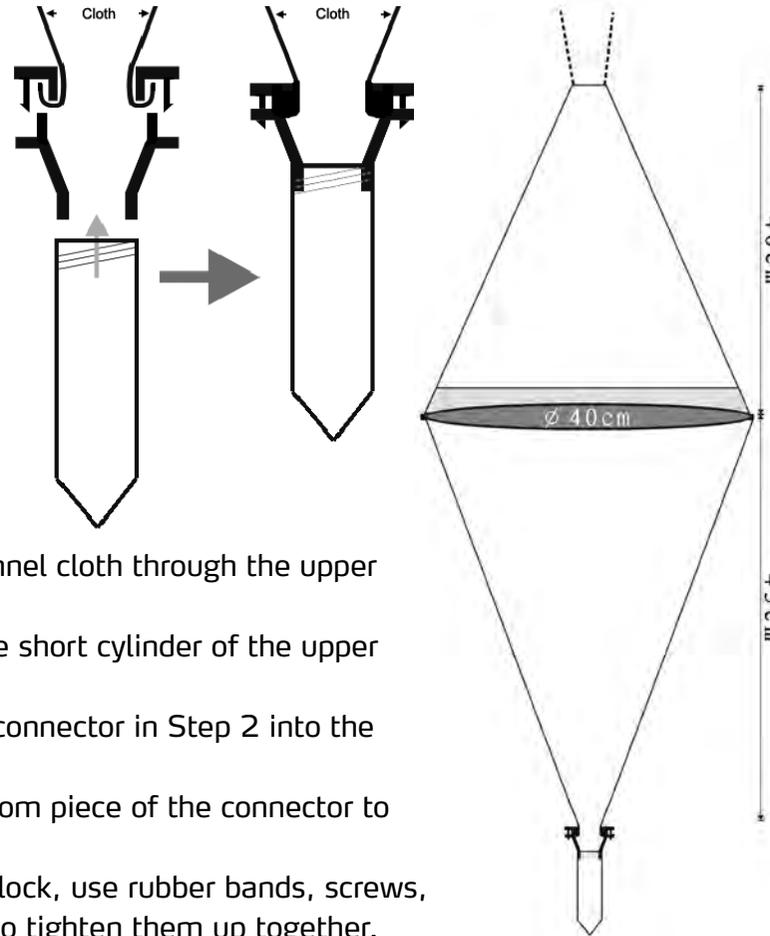

Pic. A

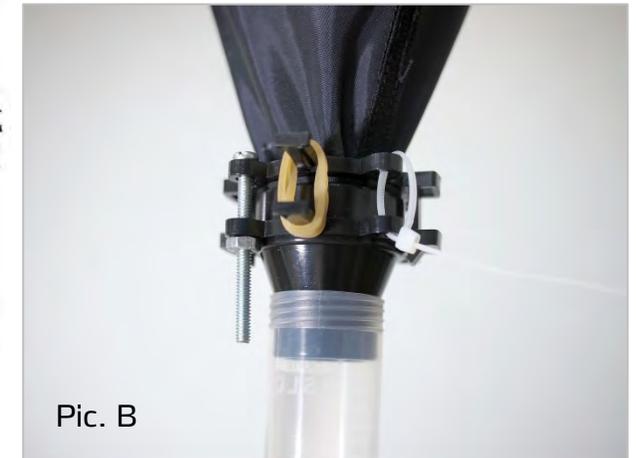

Pic. B

Have any questions? Please email us: [bugdorms@megaview.com.tw](mailto:bugdorms@megaview.com.tw) or [bugdorms@gmail.com](mailto:bugdorms@gmail.com)